

Product Name	Cat. No.	Pack Size
--------------	----------	-----------


Pre-Stained Protein Ladder	# BB-PS20	250 µl
----------------------------	-----------	--------

Pre-Stained Protein Ladder (Cat # BB-PS20) comes in a ready to use loading formulation. Pack size 250 µl (50 Loadings or more).

Maximum loading volume of this ready to load Pre-Stained Protein Ladder is 5 µl in a standard PAGE.

**Description:**

BioBharati Pre-stained Protein Ladder is a mixture of seven highly purified recombinant proteins covalently coupled to a dye that resolves into 7 sharp bands when electrophoresed. The protein concentrations are carefully balanced for even intensity. The orange 35 KDa band serve as the reference indicators. The apparent molecular weights of the Pre-stained Protein Ladder bands were determined on Tris-glycine SDS-PAGE gels by comparison to the unstained Protein Ladder. This Pre-stained Protein Ladder is suitable to use in SDS-PAGE and western blotting applications. It allows continuous monitoring of protein separations during electrophoresis and also provides a quick and easy way to assess blotting efficiency.

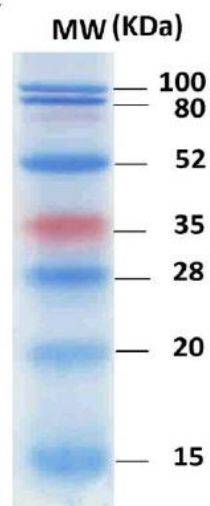

Pre-Stain Protein Ladder (Cat # BB-PS20) in a Western Blot after 12.5%, Tris-Glycine SDS-PAGE.

**Storage Condition: Store at -20°C.**

The Pre-stained Protein marker comes ready for direct loading on a SDS-PAGE. No need to do any other processing for loading in a SDSPAGE. Mix well (vortex) and directly load on the gel.

The Pre-stained Protein marker should be aliquoted into multiple tubes to prevent repeated freeze & thaw.