

Product Name	Cat. No.	Pack Size
Protein Ladder	Cat # BB-P10	500 µl
	Cat # BB-P11	1000 µl

Product type: Protein Molecular Weight ladder.

Description: Protein molecular weight marker is composed of eight recombinant proteins. Protein Concentrations for each protein are 0.1 - 0.2 mg/ml. This protein standard is ready to load in SDS-PAGE and give eight distinct bands on staining with Coomassie Brilliant Blue.


Note: Don't freeze thaw the whole Pack repeatedly. Mix well before use.

Loading: Mini Gel applications 5 µl / well

Large Gel applications 10 µl / well

Storage buffer: 25 mM Tris-HCl Buffer pH 7.4, containing 1 mM EDTA and 40 mM DTT.

Storage instructions: Store at -20°C (Recommended).

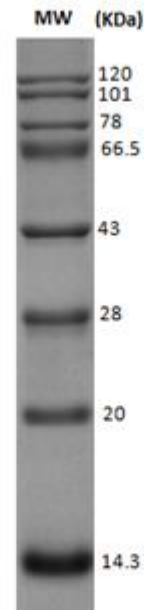

12.5% Tris-glycine SDS-PAGE Staining with Coomassie Brilliant Blue